



TECHNICAL PROTOCOL
FOR
704-Cre; Amp^R
expression plasmid
(A114)

Gene Bridges GmbH
Im Neuenheimer Feld 584
69120 Heidelberg, Germany
Tel + 49 (0)6221 1370811
Fax + 49 (0)6221 1370829
Email: products@genebridges.com
www.genebridges.com

CONTENTS

1 Eppendorf tubes + manual

1. 704-Cre; Amp^R: expression plasmid for Cre recombinase
(0.2 µg/µl, 20 µl)
2. This manual

Store tube at -20°C

Please read

The products listed in this manual are for research purposes only. They are not designed for diagnostic or therapeutic use in humans, animals or plants. The Red[®]/ET[®] recombination technology is the intellectual property of Gene Bridges GmbH.

Conditions of use

3.1 Purchaser will not manufacture, copy, reproduce, transmit, distribute, sell, lease, transfer, or improve upon the MATERIALS without prior written consent from GENE BRIDGES.

3.2 All MATERIALS relating Technologies shall be purchased from GENE BRIDGES or its authorized distributors. Use of any of the stated products from a source other than GENE BRIDGES will exempt GENE BRIDGES from any and all liabilities and warranties.

3.3 All MATERIALS purchased by research organizations, universities and other non-profit organizations may not be used for any commercial purpose. These MATERIALS are to be used for research purposes only. The MATERIALS may not be used to provide a commercial or non-commercial service, of any kind.

3.4 A purchase of MATERIALS by a private consumer is neither intended nor permitted.

Short Description:

704-Cre; Amp^R plasmid is designed for use in Cre-mediated genomic manipulations. The plasmid has a pSC101 origin which maintains low copy and replicates at 30°C. The plasmids will not propagate and will get lost when incubated at 37°C.

The expression of the Cre-recombinase is driven by the thermosensitive promoter cI578 (λ_{PR} promoter). Therefore, the expression of Cre is repressed at 30°C and induced between 37-42°C.

The plasmid carries an ampicillin resistance.

Note:

704-Cre; Amp^R plasmid was sequenced completely.

Reference:

Buchholz, F., Angrand, P.-O. and Stewart, A.F. (1996) „A simple assay to determine the functionality of Cre or FLP recombination targets in genomic manipulation constructs” *Nucleic Acids Research* 24, 3118-3119.

Zhang, Y., Buchholz, F., Muyrers, J.P.P. and Stewart, A.F. (1998) “A new logic for DNA engineering using recombination in *Escherichia coli*” *Nature Genetics* 20, 123-128.


Site Specific Recombination to Remove Selection Marker.

1. 704-Cre; Amp^R plasmid is transformed into an *E.coli* strain, which contains a targeting plasmid carrying a floxed selection marker (e.g. pPGK-neo resistance gene).
2. After transformation (electroporation or heat shock), add 1 ml of LB medium to the tube and incubate at 30°C for 1.5 h with shaking.
3. Streak out the cells on LB plates containing 50 µg/ml of ampicillin (amp) plus selection marker for the targeting plasmid.
4. Incubate at 30°C for more than 24 hours (since the colonies grow slowly).
5. Pick a single colony and grow the cells in 1 mL of LB medium containing the antibiotic of the targeting plasmid in a suitable concentration at 30°C for 2-3 hours.
6. Incubate over night at 37°C

(During incubation at 37°C, Cre protein is expressed and the loxP sites recombined; at the same time 704-Cre; Amp^R plasmid is lost.)
7. Prepare plasmid DNA and digest part of the DNA to check the restriction pattern.
8. Re-transform the checked DNA to remove of the un-recombined plasmid.


About 95% of the floxed fragment will be recombined. Step 8 is therefore important to obtain the pure and recombined plasmid.

Map:

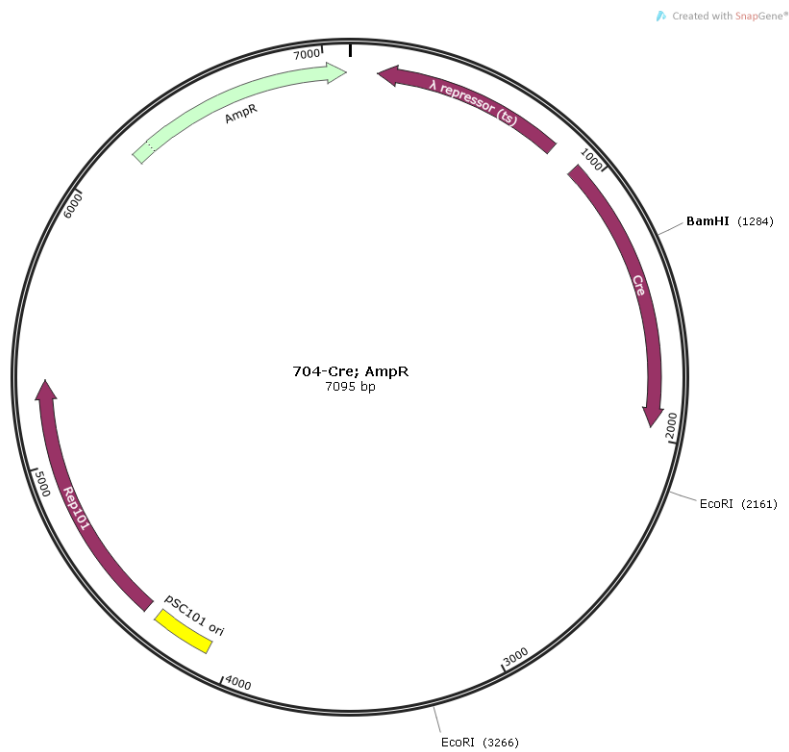

Restriction pattern of plasmid 704-Cre; Amp^R

<i>Bam</i> HI digestion:	7095 bp
<i>Eco</i> RI digestion:	1105 bp and 5990 bp
<i>Bam</i> HI and <i>Eco</i> RI digestion:	877 bp, 1105 bp and 5113

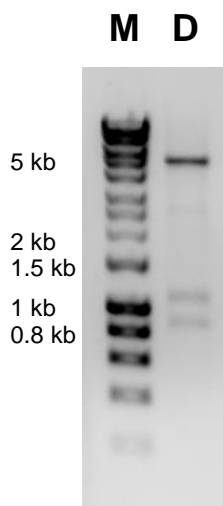

M: HyperLadder™ 1 kb, Bioline
 D: Digestion of 704-Cre; Amp^R with *Bam*HI and *Eco*RI