



TECHNICAL PROTOCOL

FOR

pCAGGS-FLPe

Enhanced Eukaryotic FLP Expression
Plasmid

(A201, A202)

CONTENTS

1 Eppendorf tubes + manual

1. pCAGGS-FLPe: expression plasmid for FLPe recombinase (0.2 µg/µl, 20 µl)
2. This manual

Store tube at -20°C

Please read

The products listed in this manual are for research purposes only. They are not designed for diagnostic or therapeutic use in humans, animals or plants. The Red[®]/ET[®] recombination technology is the intellectual property of Gene Bridges GmbH.

Conditions of use

3.1 Purchaser will not manufacture, copy, reproduce, transmit, distribute, sell, lease, transfer, or improve upon the MATERIALS without prior written consent from GENE BRIDGES.

3.2 All MATERIALS relating Technologies shall be purchased from GENE BRIDGES or its authorized distributors. Use of any of the stated products from a source other than GENE BRIDGES will exempt GENE BRIDGES from any and all liabilities and warranties.

3.3 All MATERIALS purchased by research organizations, universities and other non-profit organizations may not be used for any commercial purpose. These MATERIALS are to be used for research purposes only. The MATERIALS may not be used to provide a commercial or non-commercial service, of any kind. The use of the CAGGS promoter is covered by patents (CST Research Institute, Kumamoto, Japan).

3.4 A purchase of MATERIALS by a private consumer is neither intended nor permitted.

Short Description:

pCAGGS-FLPe-IRESpuro is an improved vector for expression of an enhanced FLP recombinase in mouse ES cells and oocytes. FLPe shows enhanced thermo stability at 37°C compared to the normal FLP recombinase (see figure 1). After micro-injection of pCAGGS-FLPe into oocytes approximately one-third of heterozygotic mice born show complete FLP recombination (Schaft J. et al., Genesis 31, 6-10 (2001)).

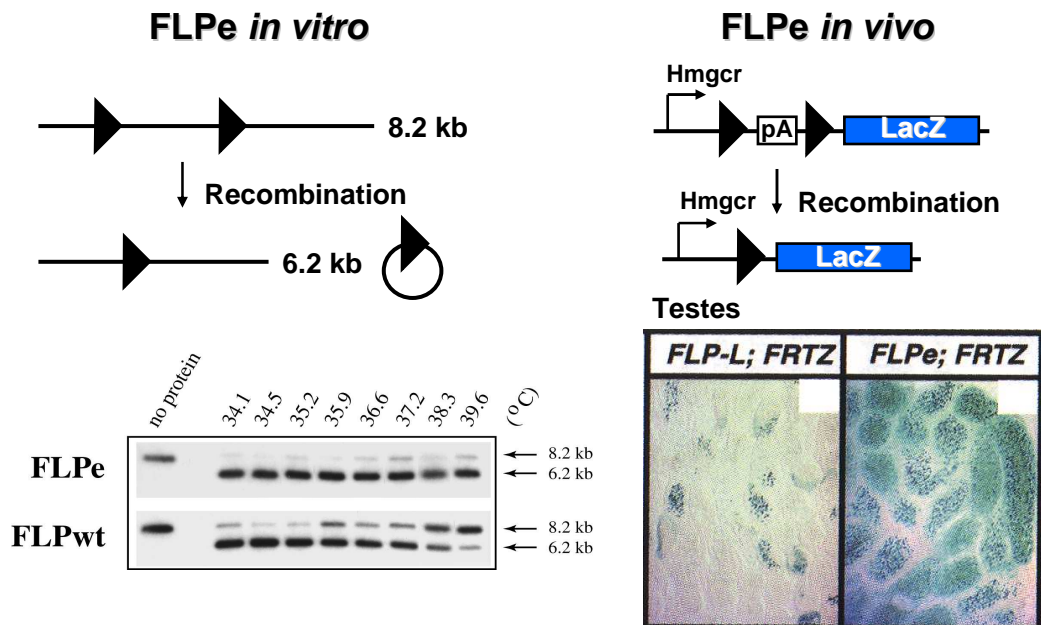


Fig. 1: Comparison of FLPe and FLPwt *in vitro* and *in vivo*. Effect of different temperatures on the recombination efficiency of FLPe in comparison to FLPwt (left part). Treatment with FLPe instead of FLPwt shows a much higher percentage of recombined products when incubated at 37°C. FLPe deleter strains achieve complete excision of the polyA signal in the germ line in comparison to FLP-L strain (for details see Rodriguez et al. 2000).

pCAGGS-FLPe-IRESpuro carries a strong, constitutive expression cassette for an *in vitro*-evolved FLP recombinase (FLPe; Buchholz F., Angrand P.O. and Stewart A.F. Nature Biotechnology 16, 657-662 (1998)) under the control of the chicken- β -actin promoter and an hCMV immediate early enhancer. The use of the chimeric CMV enhancer/ β -actin promoter leads to a ubiquitous expression profile in eukaryotes. The FLPe recombinase carries a minimal, nuclear localization signal from the SV40 large T antigen and is operatively linked by an IRES to a puromycin resistance gene. Additional elements of the expression cassette are an intron with splice donor and acceptor sites upstream of the FLPe gene and a bovine growth hormone polyadenylation signal downstream of the puromycin resistance gene.

The pCAGGS-FLPe plasmid is appropriate for efficient excision of DNA stretches flanked by FRT sites e.g. for excision of a resistance cassette in a conditional allele in eukaryotic cells. The plasmid carries an ampicillin resistance cassette.

Electroporation in ES cells:

10 μ g of supercoiled pCAGGS-FLPe plasmid were used for a single electroporation of 5×10^6 cells in 500 μ l. Cells were diluted 1000-fold afterwards and aliquots plated without selection. 22 hours later cells were changed to 1 μ g/ml puromycin. Cells were left under puromycin selection for 48 hours and changed afterwards to media without puromycin. (see also: Schaft et al. 2001). Using transient puromycin selection, some 10% of all colonies show recombination (without replating or mosaics).

Literature:

- Brown E.J. and Baltimore D. 2003: Essential and dispensable roles of ATR in cell cycle arrest and genome maintenance. *Genes & Development* 17: 615-628.
- Buchholz F., Angrand P.O. and Stewart A.F. 1998: Improved properties of FLP recombinase evolved by cycling mutagenesis. *Nature Biotechnology* 16: 657-662.
- Cope D.W. et al. 2004: Abolition of zolpidem sensitivity in mice with a point mutation in the GABA_A receptor γ 2 subunit. *Neuropharmacology* 47: 17-34.
- Morey C. et al. 2004: The region 3' to Xist mediates X chromosome counting and H3 Lys-4 dimethylation within the Xist gene. *The EMBO J.* 23: 594-604.
- Rodriguez C. et al. 2000: High-efficiency deleter mice show that FLPe is an alternative to Cre-*loxP*. *Nature Genetics* 25: 139-140.
- Schaft J. et al. 2001: Efficient FLP Recombination in mouse ES cells and oocytes. *Genesis* 31: 6-10.
- Umana P. et al. 2001: Efficient FLPe recombinase enables scalable production of helper-dependent adenoviral vectors with negligible helper-virus contamination. *Nature Biotechnology* 19: 582-585.
- Yadav et al. 2003: Specific protein methylation defects and gene expression perturbations in coactivator-associated arginine methyltransferase 1-deficient mice. *PNAS* 100: 6464-6468.

Map:

